DE

EN

FR

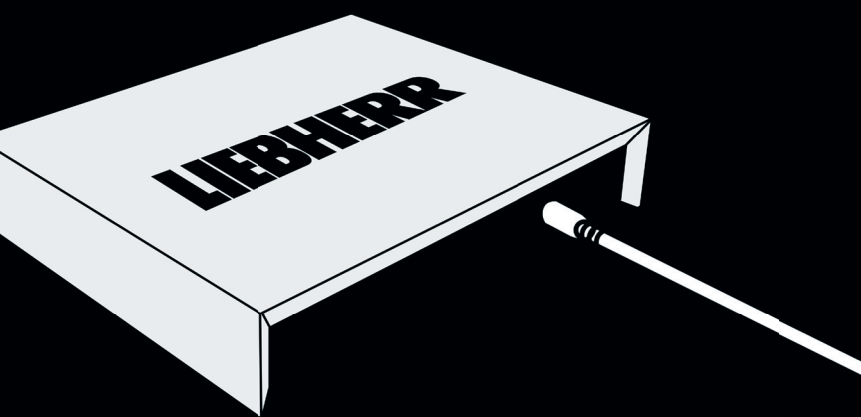
ES

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Quick Start Guide

Guide rapide

Guía rápida

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- For assistance with set-up, please contact Liebherr Service (see contact info).
- The plug-in power supply must be plugged together for the local socket standard. Select an adapter and slide onto the power supply unit until it clicks into place. A power connection designed for the power supply unit should be located near the SmartCoolingHub.
- Broadband Internet connection with wireless network (WLAN 2.4 GHz) or wired (Ethernet/LAN)
- The SSID must be visible for the WLAN connection.
- Share the following network ports within your network and at the Internet transition point (keyword: Firewall):
 - DNS: Port 53(OUT)UDP/TCP
 - NTP: Port 123 (OUT)/UDP
 - HTTPS: Port 443 (OUT)/TCP
 - Secure MQTT: Port 8883 (OUT)/TCP, 8884 (OUT)/TCP

WiFi IEEE 802.11 b/g/n 2.4 GHz, WPA2-PSK

Alternatively: (automatic) UPnP port sharing (Universal Plug and Play)

- DHCP mode active (automatic IP address assignment)
- The MAC address of the SmartCoolingHub must be shared in the network (see label on the back of the SmartCoolingHub if a MAC filter is used in your network).
- An Internet-enabled device with a current browser version and a minimum resolution of 1024x768 pixels is required to access SmartMonitoring.
- For your convenience, make a note of the serial numbers of your refrigerators in advance.
- A MyLiebherr customer profile is required to log in to the SmartMonitoring dashboard (registration possible during set-up).
- You will need additional accessories to set-up other appliances.
- Recommendation: one SmartCoolingHub per room
- If the distance between the SmartCoolingHub and the appliance is greater than 50 m, you should consult a qualified technician for installation.

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 Start the set-up at the following address:

smartmonitoring.liebherr.com/setup or scan this QR-Code:

A square QR code with a black and white pixelated pattern, used for quick access to the setup page.

You will also find the following there:

- Set-up videos
- Appliance instruction manual
- SmartMonitoring instruction manual

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IT Guida rapida

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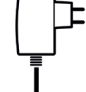
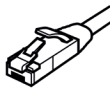
Version 1.8

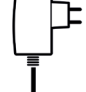
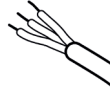
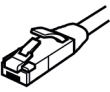
latest version available at home.liebherr.com/smartmonitoring

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